

baxterTM

research robot with SDK pre-delivery guide

This guide contains the information you need to:

- Prepare the Workspace
- Set Up the SDK
- Move and Set Up Baxter

Baxter is on its way!

We are delighted that you have chosen the Baxter Research Robot with Software Development Kit (SDK) to join your team. This document is designed to help ensure a smooth installation of Baxter in your facility, including helping you determine the optimal location and tools/equipment required for set up.

The key steps to setting up Baxter are:

Prepare the Workspace

(if mounting Baxter to a table, drill holes)

Set Up the SDK

Move Boxes

Have on hand: Pallet jack

1

Assemble the Pedestal

(if purchased)

Have on hand: 27 mm wrench, tape measure

2

Place Baxter

Mount to a pedestal or table/bench, POWER ON

Have on hand: 17 mm wrench, 1/2" wrench, tape measure, flat-head screwdriver, hoist/lift

3

Attach Grippers

Prepare the Workspace

To prepare for the arrival of Baxter:

1. Identify members of your staff to unpack and set up Baxter.
2. Ensure the unloading area and truck height are compatible, or that the truck is lift-capable.
3. Ensure any elevators or ramps are available and in place to move Baxter to its new location. Weight and dimensions of Baxter and optional pedestal, including packaging:
 - Baxter: 41 x 32 x 49 in (102 x 82 x 110 cm), 244 lb
 - Pedestal: 49 x 36 x 19 in (124 x 92 x 48 cm), 214 lb
4. Baxter is designed for easy set-up by following the simple directions included in the packaging. If you have any questions, please contact Rethink Support.
5. Baxter weighs 165 lb. If you are installing Baxter on a bench/table, identify a solid bench or table with a flat, level surface, and four sturdy legs. Anchor the bench/table securely to the floor.

Contact Rethink Robotics Customer Support at rsdk.support@rethinkrobotics.com for a copy of the bolt pattern template and mounting requirements.

6. Set up the Baxter workspace. Consider the following requirements:
 - Power: Standard 120V AC, 6A
 - Vacuum gripper (if ordered): Oil-free, compressed air @ 60 PSI min to 90 PSI max
 - Workspace dimensions: see “Baxter dimensions” on pages 5 through 8 to ensure you allow sufficient workspace to accommodate Baxter.

Set Up the SDK

If you do not already have a github account please create one at github.com.

Please email Research SDK support at rsdk.support@rethinkrobotics.com with your github username, PO/order number, and serial number of your robot. We will get back to you shortly with full access to the Wiki, which will explain how to set up the Research SDK with your Baxter.

Move Boxes and Set Up Baxter

When the shipment arrives:

1. Before signing the delivery sheet, confirm all parts ordered are included in the shipment, and all packaging is intact.
2. Prepare members of your staff to unpack and set up Baxter. Set-up instructions are located in the accessories box. You will need the following tools and equipment:
 - Pallet jack
 - Hoist/lift (with chain) capable of lifting 165 lb
 - ½-in wrench
 - 17 mm wrench
 - 27 mm wrench (pedestal only)
 - Scissors
 - Flat-head screwdriver

Additional tools are provided in the Baxter and gripper shipment.

3. Move the boxes by pallet jack into the workspace area, and unpack there; or if you purchased a pedestal, and the path from the delivery area to the Baxter workspace is flat (no thresholds, etc.), you may prefer to unpack Baxter at the delivery area, mount it onto the pedestal, and roll it to the work location.


Note that the pedestal weighs 134 lb, and Baxter weighs 165 lb. Follow proper lift guidelines and restrictions for these weights.

4. If attaching Baxter to a prepared bench or table, use four M10 bolts of sufficient length.
5. Retain all packing material for any future service needs.


Unpacking and set-up should take under an hour.

Baxter dimensions



Top view

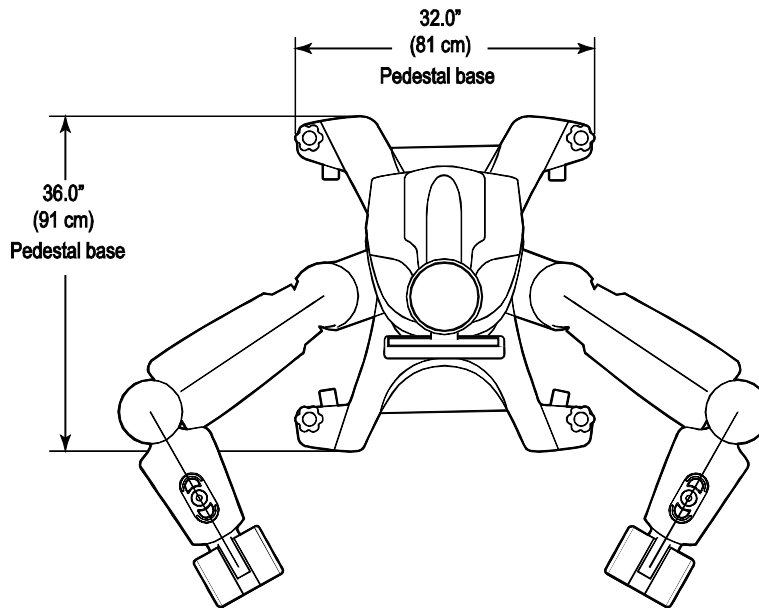


Front view

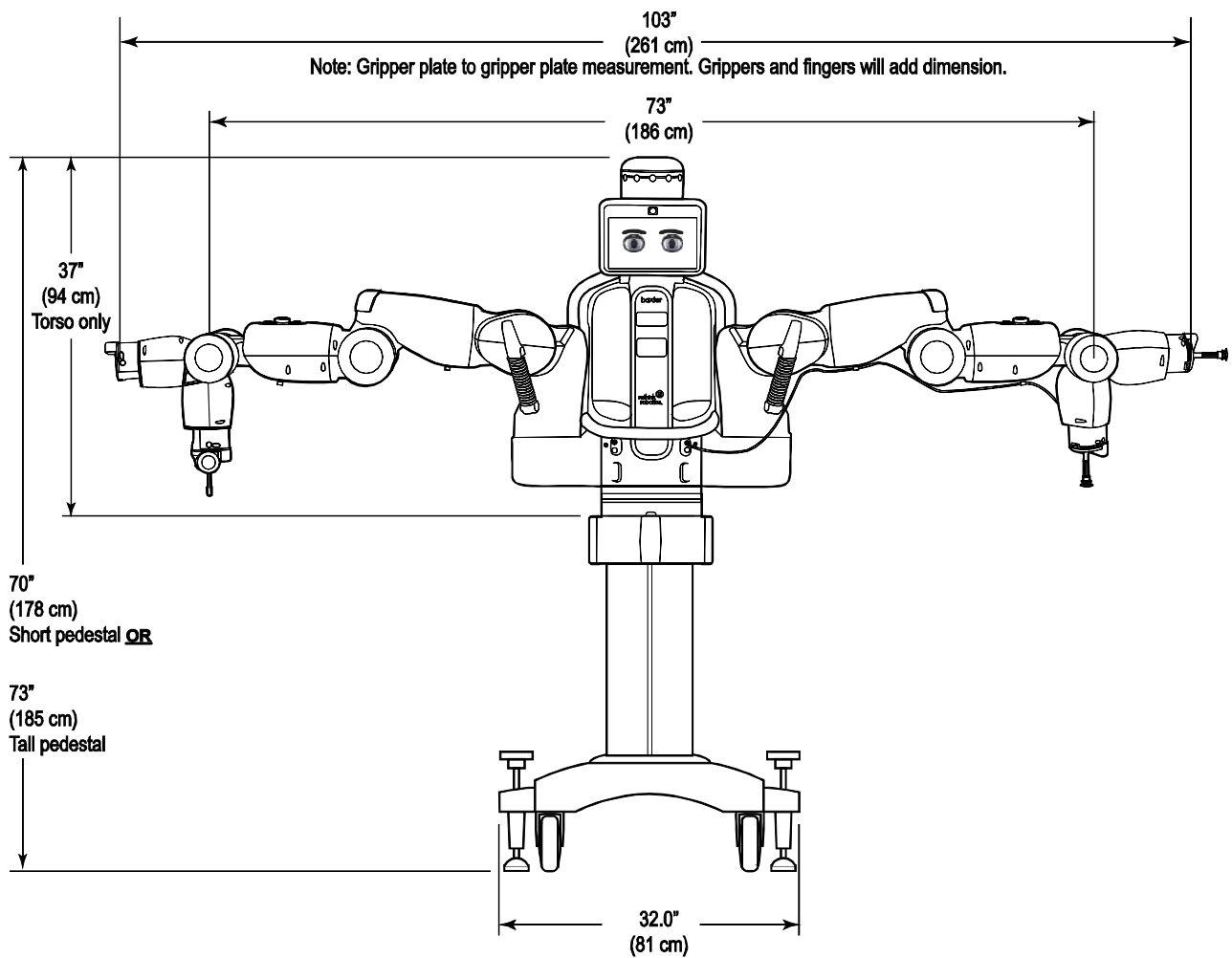



Side view

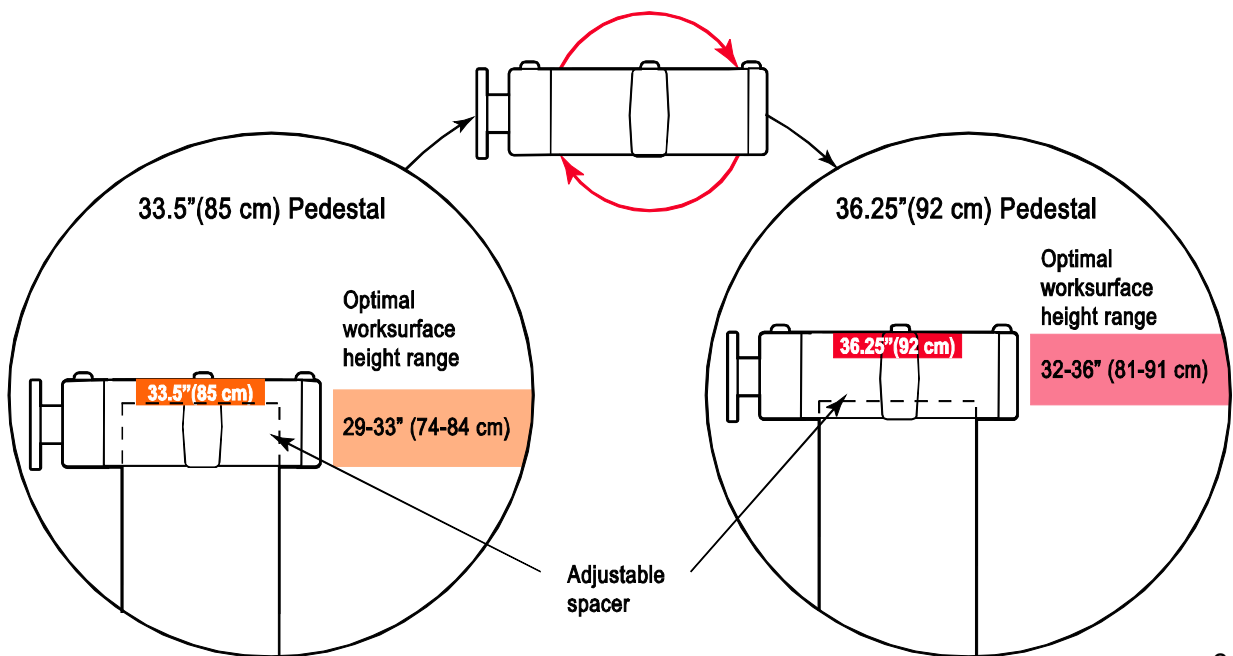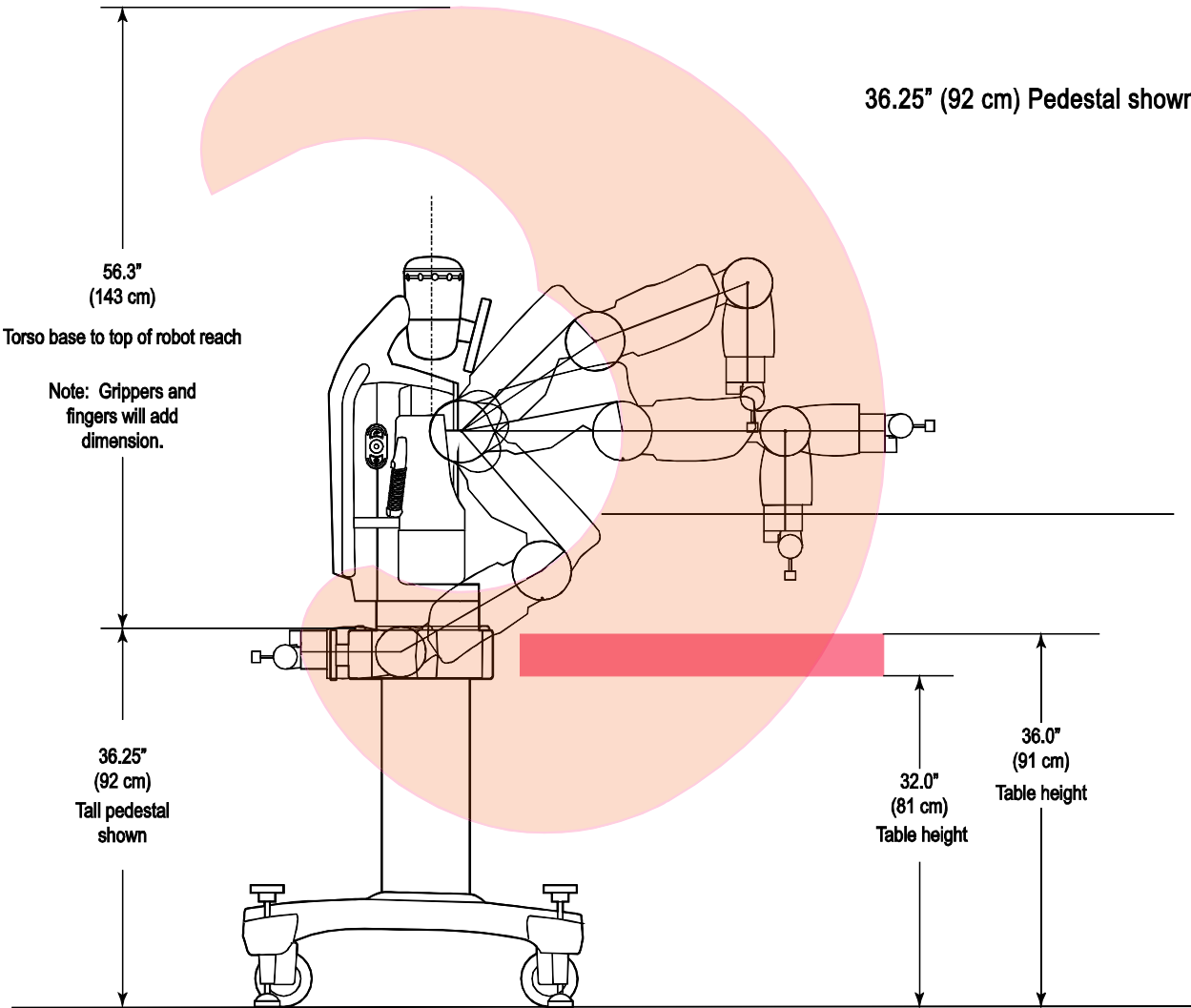

Top view, arms extended

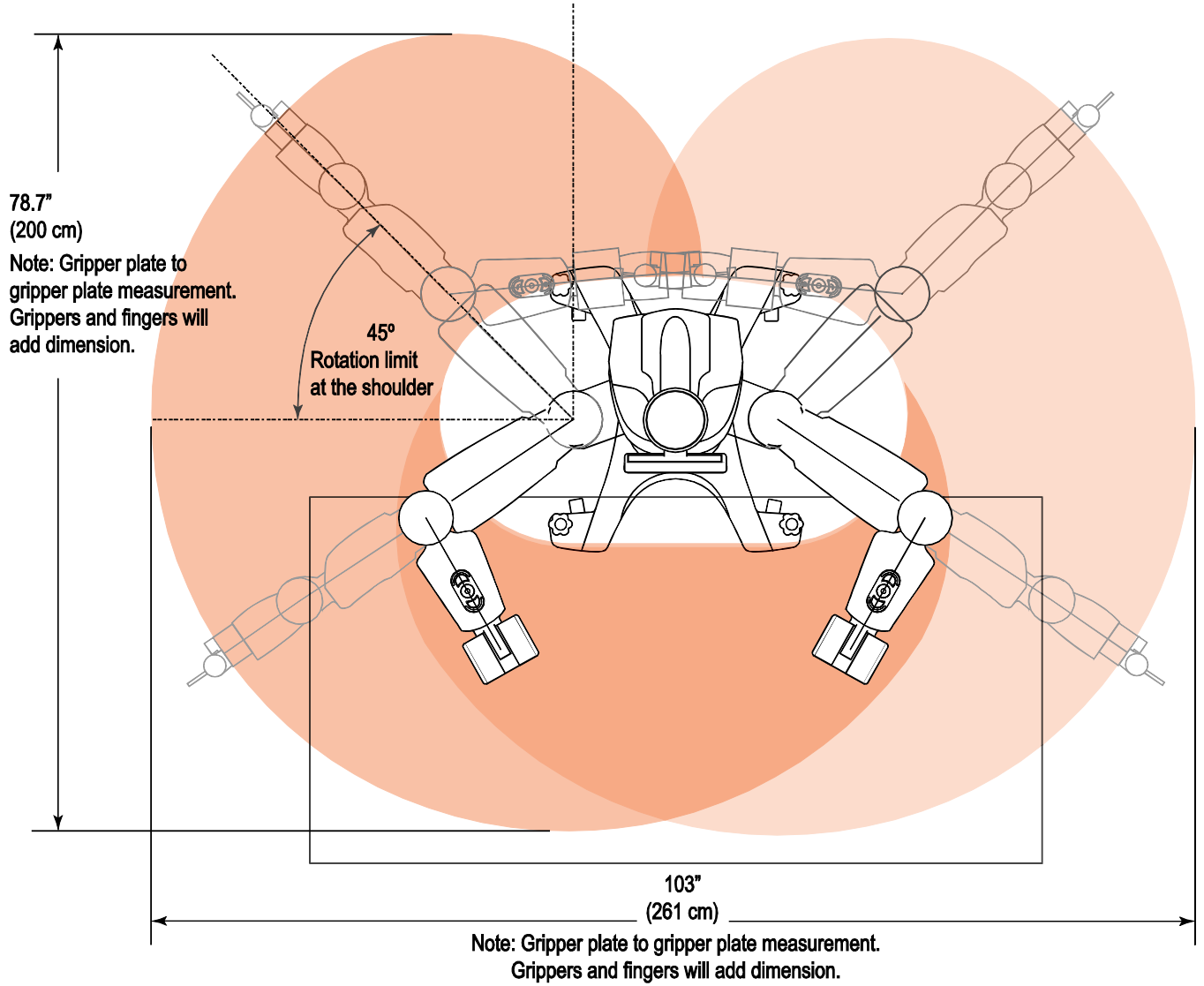

A bit about Baxter

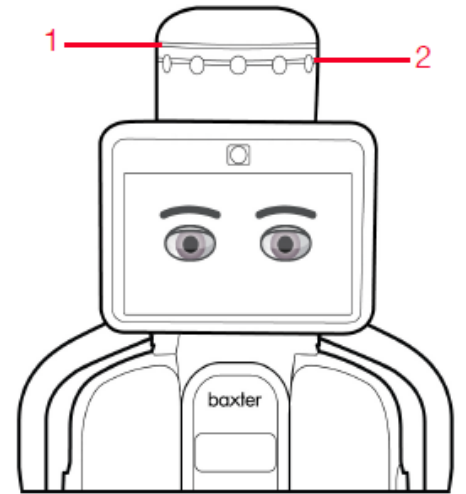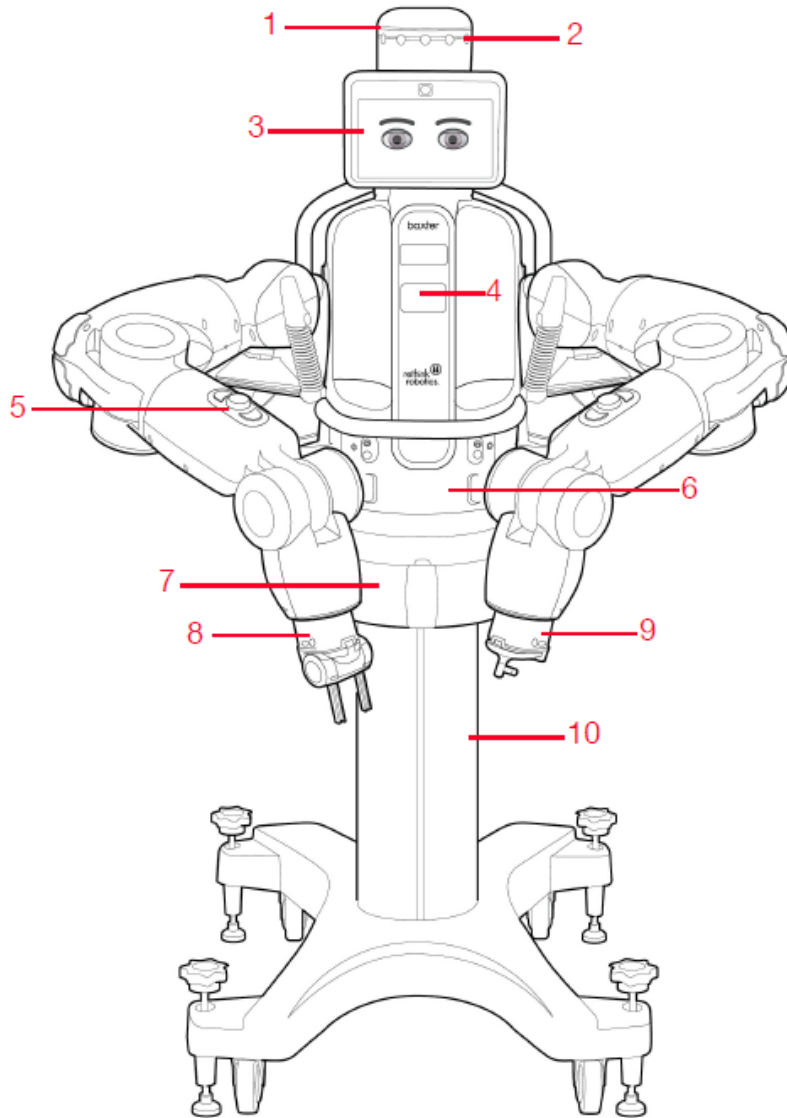

1. Condition ring
2. Attention ring
3. Display
4. Work lights
5. Navigator
6. Lower front panel
7. Waist
8. Training cuff with parallel gripper
9. Training cuff with vacuum gripper
10. Pedestal



©2013 Rethink Robotics, Inc. 27-43 Wormwood St, Boston, MA 02210. All Rights Reserved.
Rethink Robotics and Baxter are trademarks of Rethink Robotics, Inc.
U.S. Patents Pending.